

# Drinking Water Sample Set 2

## LEAD AND COPPER

Collect sample after water has sat motionless in pipes for at least 6 hours!

## DBP's

Disinfection By-Products  
No air in vials once they are filled.  
Do Not Rinse!

Gross Alpha, Uranium,  
and Combined:  
Radium 226 & Radium 228



Trip Blank  
(Do not open)



Add 5 drops from the blue capped dropper to each of the 524 vials.

Preservative already in the 552 vials.

Add both red capped vials to each Container

V.4/11

Wyoming

Always fill the sample container first, then add the preservative

[www.energylab.com](http://www.energylab.com)

