

# Drinking Water Sample Set 1

**IOCs - Inorganic Compounds**  
Metals    Cyanide    Fluoride

**VOCs 524.2**  
Volatile Organic Compounds  
No air in vials once they are filled.  
Do Not Rinse! (Add 5 drops from the blue capped dropper to each vial)

Add contents of red capped vial.

Add contents of green capped vial.

**Trip Blank**  
(Do Not Open)

No air in vials once they are filled. Do Not Rinse!

**SOCs - Synthetic Organic Compounds**

504	547	548.1	549.2	515.4	525.2	531.1
EDP, DBCP Pesticides	Herbicides	Semi-Volatile Compounds	Pesticides	Herbicides	Semi-Volatile Compounds	Carbomate Pesticides

Note: All of the above SOC bottles are pre-preserved, DO NOT RINSE!



No air in vials once they are filled. Do Not Rinse!

Add contents of blue capped vial to each bottle.

V. 12/14